

AssayMax™ Human Lactoferrin Autoantibody ELISA Kit (anti-Lactoferrin IgG)

Assaypro LLC 3400 Harry S Truman Blvd St. Charles, MO 63301 T (636) 447-9175 F (636) 395-7419 www.assaypro.com

For any questions regarding troubleshooting or performing the assay, please contact our support team at support@assaypro.com.

Thank you for choosing Assaypro.

Assay Summary

Step 1. Add 50 μ l of Standard or Sample per well. Incubate 2 hours.

Step 2. Wash, then add 50 μl of Biotinylated Antibody per well. Incubate 1 hour.

Step 3. Wash, then add 50 μl of SP Conjugate per well. Incubate 30 minutes.

Step 4. Wash, then add 50 μ l of Chromogen Substrate per well. Incubate 25 minutes.

Step 5. Add 50 μ l of Stop Solution per well. Read at 450 nm immediately.

Symbol Key

Consult instructions for use.

Assay Template

12								
11								
10								
6								
∞								
7								
9								
ß								
4								
æ								
2								
1								
	A	B	C	D	Е	Ŀ	Ð	т

AssayMax[™] Human Lactoferrin Autoantibody ELISA Kit (anti-Lactoferrin IgG)

Catalog No. EL7011-1 Sample insert for reference use only

Introduction

Lactoferrin is an 80 kDa iron-binding glycoprotein produced by many exocrine glands with a major constituent in the secondary granules of neutrophilic leukocytes. Serum lactoferrin concentration is much higher during inflammation (1). Lactoferrin is known to be an immune modulator or enhancer due to specific receptors for lactoferrin that are found on many key immune cells, such as lymphocytes, monocytes, and macrophages. Lactoferrin is known to be directly involved in the up-regulation of natural killer (NK) cell activity (2). Lactoferrin is present in maternal milk, saliva, tears, vaginal secretions, semen, bronchoalveolar lavage fluid, and specific granules of polymorphonuclear leukocytes (PMNs) (3). Lactoferrin is found mainly in the oral cavity where it can come into direct contact with pathogens, such as viruses, bacteria, etc. Lactoferrin directly inhibits viruses by binding to viral receptor sites, thus preventing the virus from infecting healthy cells. Lactoferrin has a direct bactericidal function to certain bacteria, such as Streptococcus mutans, Vibrio cholerae, Escherichia coli, Actinobacillus actinomycetemcomitans, and Legionella pneumophila (2-4). Also, it has a bacteriostatic effect that deprives iron-requiring bacteria of this essential growth nutrient (4). Lactoferrin is also considered an antioxidant that scavenges free iron, helping to prevent uncontrolled iron based free radical reactions, thus protecting certain cells from peroxidation (2). Autoantibodies against lactoferrin belong to the pANCA class (Perinuclear Anti-Neutrophil Cytoplasmic Antibodies). Anti-lactoferrin antibodies are found in patients with rheumatoid arthritis with vasculitis (5) as well as in reactive arthritis and ankylosing spondylitis (6).

Principle of the Assay

The AssayMax[™] Human Lactoferrin Autoantibody ELISA (Enzyme-Linked Immunosorbent Assay) Kit is designed for the quantitative determination of autoimmune response (IgG) to a target antigen (lactoferrin). The kit detects autoantibodies in **plasma and serum samples**. This assay employs a quantitative **sandwich enzyme immunoassay technique that measures autoantibodies (anti-Lactoferrin IgG) in approximately 4 hours**. A lactoferrin antigen has been pre-coated onto a 96-well microplate with removable strips. An autoantibody specific for lactoferrin in standards and samples is sandwiched by the immobilized antigen and a biotinylated polyclonal antibody specific for human IgG, which is recognized by a streptavidin-peroxidase (SP) conjugate. All unbound material is washed away and a peroxidase enzyme substrate is added. The color development is stopped and the intensity of the color is measured.

Caution and Warning

- This product is for **Research Use Only** and is not intended for use in diagnostic procedures.
- Prepare all reagents (diluent buffer, wash buffer, standard, biotinylated antibody, and SP conjugate), as instructed, prior to running the assay.
- Prepare all samples prior to running the assay. The dilution factors for the samples are suggested in this insert. However, the user should determine the optimal dilution factor.
- Spin down the SP conjugate vial and the biotinylated antibody vial before opening and using contents.
- The Stop Solution is an acidic solution.
- The kit should not be used beyond the expiration date.

Reagents

- Human Lactoferrin Autoantibody Microplate: A 96-well polystyrene microplate (12 strips of 8 wells) coated with a lactoferrin antigen.
- Sealing Tapes: Each kit contains 3 precut, pressure sensitive sealing tapes that can be cut to fit the format of the individual assay.
- Human Lactoferrin Autoantibody Standard: Plasma standard (120 AU, lyophilized).
- Autoimmune Biotinylated Human IgG Antibody (40x): A 40-fold concentrated biotinylated polyclonal antibody against human IgG (150 μl).
- MIX Diluent Concentrate (10x): A 10-fold concentrated buffered protein base (30 ml).
- Wash Buffer Concentrate (20x): A 20-fold concentrated buffered surfactant (30 ml, 2 bottles).
- SP Conjugate (100x): A 100-fold concentrate (80 µl).
- **Chromogen Substrate (1x)**: A stabilized peroxidase chromogen substrate tetramethylbenzidine (7 ml).
- Stop Solution (1x): A 0.5 N hydrochloric acid solution to stop the chromogen substrate reaction (11 ml).

Storage Condition

- Upon arrival, immediately store components of the kit at recommended temperatures up to the expiration date.
- Store Microplate, SP Conjugate, and Biotinylated Antibody at -20°C.
- Store Diluent Concentrate (10x), Wash Buffer, Stop Solution, and Chromogen Substrate at 2-8°C.
- Unused microplate wells may be returned to the foil pouch with the desiccant packs and resealed for up to 30 days at -20°C.
- Store Standard at 2-8°C before reconstituting with Diluent and at -20°C after reconstituting with Diluent.

Other Supplies Required

- Microplate reader capable of measuring absorbance at 450 nm
- Pipettes (1-20 µl, 20-200 µl, 200-1000 µl, and multiple channel)
- Deionized or distilled reagent grade water

Sample Collection, Preparation, and Storage

- **Plasma:** Collect plasma using one-tenth volume of 0.1 M sodium citrate as an anticoagulant. Centrifuge samples at 3000 x g for 10 minutes and collect plasma. A 40-fold sample dilution is suggested into MIX Diluent; however, user should determine optimal dilution factor depending on application needs. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles (EDTA or Heparin can also be used as an anticoagulant).
- Serum: Samples should be collected into a serum separator tube. After clot formation, centrifuge samples at 3000 x g for 10 minutes and remove serum. A 40-fold sample dilution is suggested into MIX Diluent; however, user should determine optimal dilution factor depending on application needs. The undiluted samples can be stored at -20°C or below for up to 3 months. Avoid repeated freeze-thaw cycles.

Refer to Dilution Guidelines for further instruction.

	Guidelines for Dilutions of 100-fold or Greater (for reference only; please follow the insert for specific dilution suggested)					
100x		10000x				
A)	4 μl sample : 396 μl buffer (100x) = 100-fold dilution	A) B)	4 μl sample : 396 μl buffer (100x) 4 μl of A : 396 μl buffer (100x) = 10000-fold dilution			
	Assuming the needed volume is less than or equal to 400 μl.		Assuming the needed volume is less than or equal to 400 μ l.			
	1000x		100000x			
A) B)	4 μl sample : 396 μl buffer (100x) 24 μl of A : 216 μl buffer (10x) = 1000-fold dilution	A) B) C)	4 μl sample : 396 μl buffer (100x) 4 μl of A : 396 μl buffer (100x) 24 μl of B : 216 μl buffer (10x) = 100000-fold dilution			
	Assuming the needed volume is less than or equal to 240 μl.		Assuming the needed volume is less than or equal to 240 μl.			

Reagent Preparation

- Freshly dilute all reagents and bring all reagents to room temperature before use.
- MIX Diluent Concentrate (10x): Dilute the MIX Diluent Concentrate 10fold with reagent grade water to produce a 1x solution. When diluting the concentrate, make sure to rinse the bottle thoroughly to extract any precipitates left in the bottle. Mix the 1x solution gently until the crystals have completely dissolved. Store for up to 30 days at 2-8°C.
- Human Lactoferrin Autoantibody Standard: Reconstitute the Human Lactoferrin Autoantibody Standard (120 AU) with 2.4 ml of MIX Diluent to generate a 50 AU/ml standard stock solution. Allow the vial to sit for 10 minutes with gentle agitation prior to making dilutions. Prepare duplicate or triplicate standard points by serially diluting from the standard stock solution (50 AU/ml) 4-fold with MIX Diluent to produce 12.5, 3.125, 0.781, 0.195, and 0.049 AU/ml solutions. MIX Diluent serves as the zero standard (0 AU/ml). Any remaining stock solution should be stored at -20°C and used within 30 days. Avoid repeated freeze-thaw cycles.

Standard Point	Dilution	[Lactoferrin] (AU/ml)
P1	1 part Standard (50 AU/ml)	50
P2	1 part P1 + 3 parts MIX Diluent	12.5
P3	1 part P2 + 3 parts MIX Diluent	3.125
P4	1 part P3 + 3 parts MIX Diluent	0.781
P5	1 part P4 + 3 parts MIX Diluent	0.195
P6	1 part P5 + 3 parts MIX Diluent	0.049
P7	MIX Diluent	0.0

- Autoimmune Biotinylated Human IgG Antibody (40x): Spin down the antibody briefly and dilute the desired amount of the antibody 40-fold with MIX Diluent to produce a 1x solution. The undiluted antibody should be stored at -20°C.
- Wash Buffer Concentrate (20x): Dilute the Wash Buffer Concentrate 20fold with reagent grade water to produce a 1x solution. When diluting the concentrate, make sure to rinse the bottle thoroughly to extract any precipitates left in the bottle. Mix the 1x solution gently until the crystals have completely dissolved.
- SP Conjugate (100x): Spin down the SP Conjugate briefly and dilute the desired amount of the conjugate 100-fold with MIX Diluent to produce a 1x solution. The undiluted conjugate should be stored at -20°C.

Assay Procedure

- Prepare all reagents, standard solutions, and samples as instructed. Bring all reagents to room temperature before use. The assay is performed at room temperature (20-25°C).
- Remove excess microplate strips from the plate frame and return them immediately to the foil pouch with desiccants inside. Reseal the pouch securely to minimize exposure to water vapor and store in a vacuum desiccator at -20°C.
- Add 50 μ l of Human Lactoferrin Autoantibody Standard or sample to each well. Gently tap plate to thoroughly coat the wells. Break any bubbles that may have formed. Cover wells with a sealing tape and incubate for 2 hours. Start the timer after the last addition.
- Wash the microplate manually or automatically using a microplate washer. Invert the plate and decant the contents; hit 4-5 times on absorbent material to completely remove the liquid. If washing manually, wash five times with 200 µl of Wash Buffer per well. Invert the plate each time and decant the contents; hit 4-5 times on absorbent material to completely remove the liquid. If using a microplate washer, wash six times with 300 µl of Wash Buffer per well; invert the plate and hit 4-5 times on absorbent material to completely remove the liquid.

- Add 50 μl of Autoimmune Biotinylated Human IgG Antibody to each well. Gently tap plate to thoroughly coat the wells. Break any bubbles that may have formed. Cover wells with a sealing tape and incubate for 1 hour.
- Wash the microplate as described above.
- Add 50 µl of SP Conjugate to each well. Gently tap plate to thoroughly coat the wells. Break any bubbles that may have formed. Cover wells with a sealing tape and incubate for 30 minutes. Turn on the microplate reader and set up the program in advance.
- Wash the microplate as described above.
- Add 50 μl of Chromogen Substrate to each well. Gently tap plate to thoroughly coat the wells. Break any bubbles that may have formed. Incubate in ambient light for 25 minutes or until the optimal blue color density develops.
- Add 50 µl of Stop Solution to each well. The color will change from blue to yellow. Gently tap plate to ensure thorough mixing. Break any bubbles that may have formed.
- Read the absorbance on a microplate reader at a wavelength of 450 nm **immediately**. If wavelength correction is available, subtract readings at 570 nm from those at 450 nm to correct optical imperfections. Otherwise, read the plate at 450 nm only. Please note that some unstable black particles may be generated at high concentration points after stopping the reaction for about 10 minutes, which will reduce the readings.

Data Analysis

- Calculate the mean value of the duplicate or triplicate readings for each standard and sample.
- To generate a standard curve, plot the graph using the standard concentrations on the x-axis and the corresponding mean 450 nm absorbance (OD) on the y-axis. The best fit line can be determined by regression analysis using log-log or four-parameter logistic curve fit.
- Determine the unknown sample concentration from the standard curve.
- Although normal samples have been diluted 40-fold, do not multiply the value by the dilution factor. Samples with elevated levels of autoantibodies can be diluted further; for example: 80x. Account for this further dilution factor when calculating the value of the sample.

Example	Dilution Factor	Multiplication Factor For Calculating Values
Serum with normal level of anti-Lactoferrin IgG	40x	1
Serum with elevated level of anti-Lactoferrin IgG	80x	2
Serum with elevated level of anti-Lactoferrin IgG	160x	4

Typical Data

• The typical data is provided for reference only. Individual laboratory means may vary from the values listed. Variations between laboratories may be caused by technique differences.

Standard Point	AU/ml	Average OD		
P1	50	1.827		
P2	12.5	1.187		
Р3	3.125	0.611		
P4	0.781	0.259		
P5	0.195	0.111		
P6	0.049	0.057		
P7	0.0	0.033		
Normal Level Sample (40x): Serum with normal level of anti-Lactoferrin IgG				
Elevated Level Serum with elevated lev	1.278			

Standard Curve

• The curve is provided for illustration only. A standard curve should be generated each time the assay is performed.

Reference Value

 Plasma and serum samples from healthy adults were tested (n=20). Moreover, patient serum samples containing high levels of anti-Lactoferrin IgG were tested (n=9). The following ranges have been established:

Sample	anti-Lactoferrin IgG (AU/ml)	
Normal Level	< 15	
Elevated Level	≥ 15	

• It is recommended that each laboratory establish its own normal and pathological ranges of antibodies.

Performance Characteristics

- The minimum detectable dose of anti-Lactoferrin IgG as calculated by 2SD from the mean of a zero standard was established to be 0.03 AU/ml.
- Intra-assay precision was determined by testing three serum samples twenty times in one assay.
- Inter-assay precision was determined by testing three serum samples in twenty assays.

	Intra-Assay Precision			Inter-Assay Precision		
Sample	1	2	3	1	2	3
n	20	20	20	20	20	20
CV (%)	4.7%	6.8%	5.3%	9.6%	11.5%	10.5%
Average CV (%)	5.6%				10.5%	

Troubleshooting

Issue	Causes	Course of Action		
	Use of improper	 Check the expiration date listed before use. 		
	components	 Do not interchange components from different lots. 		
		 Check that the correct wash buffer is being used. 		
		 Check that all wells are empty after aspiration. 		
	Improper wash step	 Check that the microplate washer is dispensing properly. 		
		 If washing by pipette, check for proper pipetting 		
Ę		technique.		
Low Precision	Splashing of reagents while loading wells	 Pipette properly in a controlled and careful manner. 		
re	Inconsistent volumes	 Pipette properly in a controlled and careful manner. 		
~ ~	loaded into wells	 Check pipette calibration. 		
6		 Check pipette for proper performance. 		
_	Insufficient mixing of	 Thoroughly agitate the lyophilized components after 		
	reagent dilutions	reconstitution.		
	5	Thoroughly mix dilutions.		
		Check the microplate pouch for proper sealing.		
	Improperly sealed	Check that the microplate pouch has no punctures.		
	microplate	Check that three desiccants are inside the microplate		
	Microplate was left	pouch prior to sealing.		
al	unattended between	 Each step of the procedure should be performed uninterrupted. 		
ign	steps	uninterrupteu.		
Unexpectedly Low or High Signal Intensity	Omission of step	• Consult the provided procedure for complete list of steps.		
ligh	Steps performed in	Consult the provided procedure for the correct order.		
r ,	incorrect order			
ly Low ol Intensity	Insufficient amount of	Check pipette calibration.		
N SU	reagents added to wells	 Check pipette for proper performance. 		
nte L	Wash step was skipped	 Consult the provided procedure for all wash steps. 		
lbs –	Improper wash buffer	 Check that the correct wash buffer is being used. 		
cte	Improper reagent	Consult reagent preparation section for the correct		
be	preparation	dilutions of all reagents.		
ex	Insufficient or	 Consult the provided procedure for correct incubation 		
n	prolonged incubation	time.		
	periods			
e,		 Sandwich ELISA: If samples generate OD values higher then the bishest standard agist (D1) dilute samples 		
Deficient Standard Curve Fit		than the highest standard point (P1), dilute samples		
Deficient ndard Cui Fit	Non-optimal sample	further and repeat the assay.Competitive ELISA: If samples generate OD values lower		
Fici Fit	dilution	than the highest standard point (P1), dilute samples		
Je Jd	anation	further and repeat the assay.		
tar		User should determine the optimal dilution factor for		
Ś		samples.		
		Sumples		

	Contamination of reagents	 A new tip must be used for each addition of different samples or reagents during the assay procedure. 		
	Contents of wells evaporate	 Verify that the sealing film is firmly in place before placing the assay in the incubator or at room temperature. 		
	Improper pipetting	 Pipette properly in a controlled and careful manner. Check pipette calibration. Check pipette for proper performance. 		
	Insufficient mixing of reagent dilutions	 Thoroughly agitate the lyophilized components after reconstitution. Thoroughly mix dilutions. 		

References

- (1) Naot D et al. (2005) Clinical Medicine & Research. 3(2):93-101.
- (2) Brink W. (October 2000) LE Magazine.
- (3) Yamauchi K et al. (1993) Infection and Immunity. 61(2):719-728.
- (4) Conneely OM. (2001) J Am Coll Nutr. 20(5 Suppl):389S-395S.
- (5) Coremans IE et al. (1993) Adv Exp Med Biol. 336:357-362.
- (6) Locht H et al. (1999) Clin Exp Immunol. 117(3):568.

Version 1.5